

SPECIFICATIONS

The VORTEX-GENIE® 2 Mixer is classified as "Installation Category 2"

Environmental: 0°C - 38°C (32°F - 100°F), 95% Humidity max.

| MODEL | POWER REQUIRED | AMPS |
|-----------|----------------|------|
| G560 | 120V | 0.65 |
| G560E | 230V | 0.5 |
| G560-100V | 100V | 1.0 |



ETL Listed
Conforms to UL Std. 3101-1
Certified to CAN/CSA C22.2



Weight: 4 Kg (8.8 lb)


Base Dimensions: (DxWxH) 165x122x165mm (6.5x4.8x6.5in)

PARTS ASSEMBLY LIST

To order parts for the VORTEX-GENIE 2 Mixer:

Contact your local distributor or visit www.scientificindustries.com. Please specify

Part No., quantity and electric voltage.

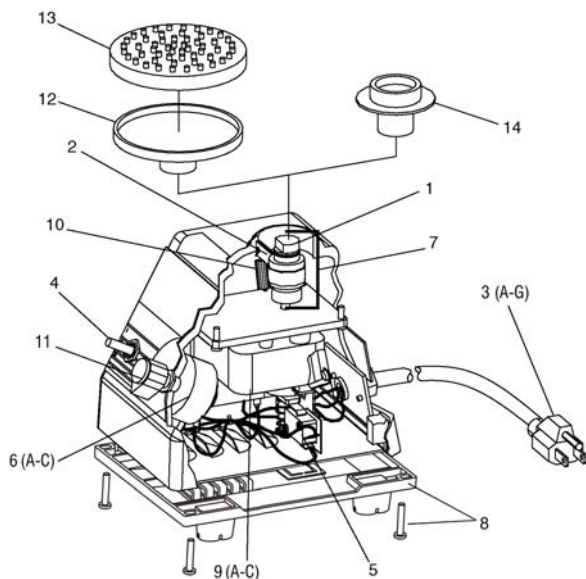

| Indicator No. | Part No. | Description | Indicator No. | Part No. | Description |
|---------------|-------------|------------------------------|---------------|-------------|----------------------------------|
| 1 | 0K-0236-902 | Bearing Retainer Kit | 6C | EB-0286-500 | Speed Control 100V |
| 2 | 0K-0236-903 | Eccentric Clamp | 7 | 0K-0236-904 | Eccentric with Clamp Assembly |
| 3A | 318-0510-02 | 120V Line Cord | 8 | 0K-0236-408 | Bottom Closure with Feet Kit |
| 3B | 0K-0246-901 | 230V Line Cord, No Plug | 9A | 0K-0236-905 | Motor 120V Assembly |
| 3C | 0K-0256-901 | 230V Line Cord, Euro Plug | 9B | 0K-0246-915 | Motor 230V Assembly |
| 3D | 0K-0266-901 | 230V Line Cord, British Plug | 9C | 0K-0286-915 | Motor 100V Assembly |
| 3E | 0K-0276-901 | 230V Line Cord, Swiss Plug | 10 | 566-0028-00 | Spring, Extension, Eccentric |
| 3F | 0K-0286-901 | 100V Line Cord | 11 | 0M-0236-209 | Knob, Push-on |
| 3G | ECP0021 | Australian Plug (only) | 12 | 0K-0500-901 | 3-inch Platform |
| 4 | ESP0004 | Switch, On/Off/On | 13 | 580-2013-00 | Rubber Cover for 3-inch Platform |
| 5 | ESP0005 | Micro Switch | 14 | 146-3011-00 | Pop-off Cup |
| 6A | EPP0005 | Speed Control 120V | | | |
| 6B | EB-0246-500 | Speed Control 230V | | | |

Vortex-Genie® is a registered trademark of Scientific Industries
© Scientific Industries Inc. 2013

U.S. Patent No. 4,781,487
065-0004-00 Rev. V

VORTEX-GENIE® 2

OPERATING INSTRUCTIONS

Models G560, G560E and G560E-100V



70 Orville Drive, Bohemia, NY 11716 U.S.A.
(631) 567-4700 • Fax: (631) 567-5896 • Toll Free: 888-850-6208
Customer Service: custsvc@scientificindustries.com
www.scientificindustries.com

VORTEX-GENIE® mixers are the "Gold Standard" for laboratory mixers. The quality and reliability, that you have come to know so well, are characteristics that are, by design, an integral part of all GENIE™ products.

APPLICATIONS

The 3-inch Platform accepts single or multiple test tubes, beakers and various flasks. The removable cover, made of thermoplastic rubber, is highly impervious to chemicals. It is, however recommended that spills are blotted quickly. Do not attempt to use the 3-inch Platform without the rubber cover.

The Pop-off Cup can accommodate tubes of virtually any size or shape from micro tubes to those up to 40mm in diameter and 200mm in length. In most instances, tubes without caps may be mixed with proper speed control. Tubes must be capped when potential aerosols may be hazardous.

Catalog numbers:

OK-0500-902 3-inch Platform with Rubber Cover
OK-0500-901 3-inch Platform without Rubber Cover
580-2013-00 Rubber Cover for 3-inch Platform

OPERATING INSTRUCTIONS

We recommend you retain the original packaging for 90 days in case you need to return the product for any reason to your distributor or Scientific Industries.

Your VORTEX-GENIE 2 Mixer is supplied with a Pop-off Cup located in the white foam box insert and a 3-inch Platform. After you unpack your VORTEX-GENIE 2 Mixer from its protective plastic bag, install either the Pop-off Cup or the 3-inch Platform by first aligning the "flat" part of the mixer shaft with the "flat" part of the attachment. **YOU MUST PRESS DOWN HARD WITH YOUR HAND UNTIL YOU HEAR THE ATTACHMENT SNAP IN PLACE.** This will ensure that the attachment will not come off the mixer during operation.

1.0 - Plug the cord into a properly grounded, three pronged outlet (**120 VAC for the G560 model, 230 VAC for the G560E model and 100V for the G560-100V model**).

2.0 - The VORTEX-GENIE® 2 Mixer has the following modes of operation:

2.1 - TOUCH MODE:

Place the toggle switch in the TOUCH position and set the speed control to any position between 3 and 10. A setting of 3 is least vigorous mixing and 10 is the most vigorous mixing. When you touch a vessel to the 3-inch Platform or Pop-off Cup, the mixer will start automatically. Releasing the pressure on the Platform or Cup will stop the mixer.



2.2 - HANDS FREE MODE:

Place the toggle switch in the ON position and set the speed control to any position between 3 and 10. Hands Free mode allows for unattended operation that will continue indefinitely until you stop the mixer by placing the toggle switch into the OFF position. In this mode, you may use any of the accessory attachments available for the VORTEX-GENIE® 2 Mixer. **Review instruction No. 3.0 before attempting to change attachments.**

Note: The speed control switch should be set between positions 3 and 10 to enable the mixing action to start. After the action has started, you may, if desired, reduce the speed below position 3 for the duration of the run. Before re-starting the mixer, re-set the speed control to a setting higher than position 3.

3.0 - To change accessory mixing attachments, the following sequence is to be followed explicitly.

3.1 - Turn the power OFF by placing the toggle switch into the OFF position.

3.2 - Grasp the installed attachment and pull straight up until it releases from the shaft. A technique of lifting the attachment with your finger tips while using your knuckles as a fulcrum will give you the leverage to remove the installed attachment component smoothly.

3.3 - Hold the part you wish to install so the "flat" part of the shaft is aligned with the "flat" of the receptacle in the attachment. Alternatively, you may place the attachment on the shaft and turn it until you feel the "flats" aligning.

3.4 - Press attachment firmly downward until it "snaps" into place. The VORTEX-GENIE 2 Mixer is now ready to use with the newly installed accessory attachment.

4.0 - With prolonged use, your VORTEX-GENIE 2 Mixer may become warm to the touch. The motor will radiate a certain amount of heat that will dissipate by the cast zinc housing.

CAUTION! Do not lift your VORTEX-GENIE 2 Mixer by its attachment. All attachments, including the Pop-off Cup and the 3-inch Platform are removable. They will pop off easily if you lift them, and snap back on again just as easily by aligning the flats and pressing down firmly.

CAUTION! Failure to follow operating instructions can compromise the user's safety.

CARE AND HANDLING

Your VORTEX-GENIE 2 Mixer should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. The finish can be washed with water (after unplugging) and soap or detergents, using a cloth or sponge. Keep the unit clean by immediately blotting any spills. Replacement parts are available through most laboratory equipment distributors or directly from Scientific Industries. Refer to "Parts Assembly List" and enclosed multi-fold brochure.

CAUTION! Unplug from power before cleaning. Do not immerse.